



## How to install an application (i.e. SpectraSuite) in a language other than the default language under which Windows XP has been installed.


1. It is assumed that any previous installation of the application has already been uninstalled.
2. Open the “Control Panel” and select “Regional and Language Options.” (If you are using the Control Panel’s “Category View,” select Date, Time, Language, and Regional Options” and then “Regional and Language Options.”)

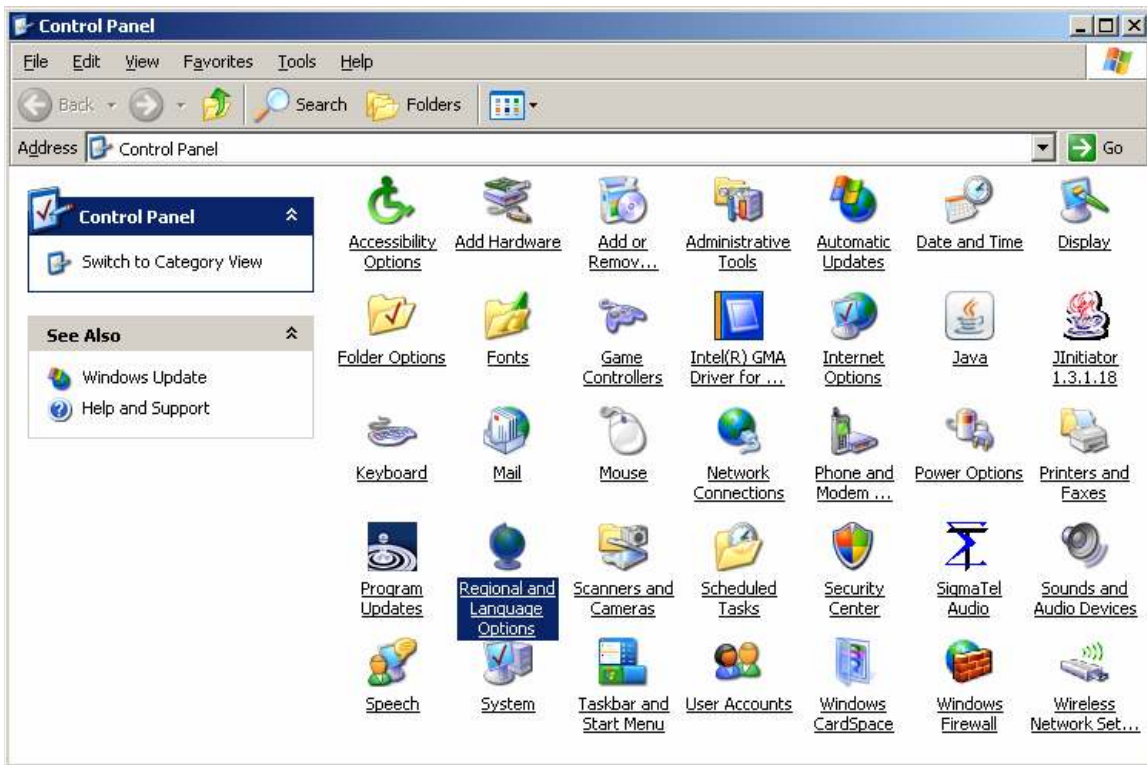

### 3. Click on the “Regional Options” tab.

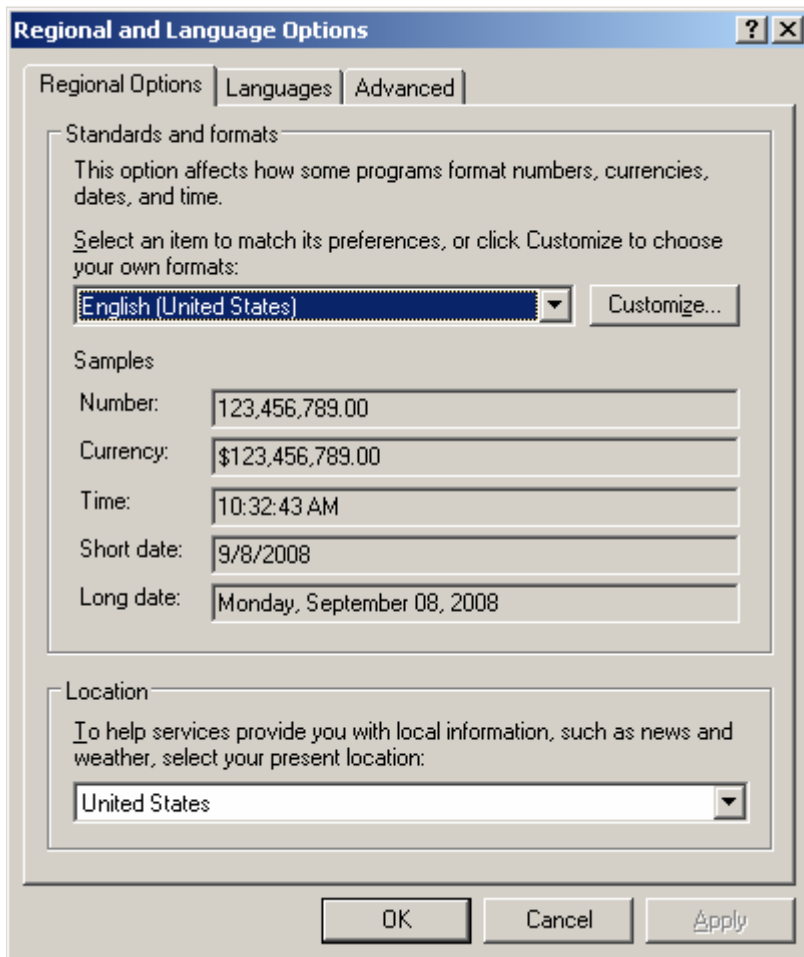

4. In the drop-down list, select the language in which you would like to install the new application.

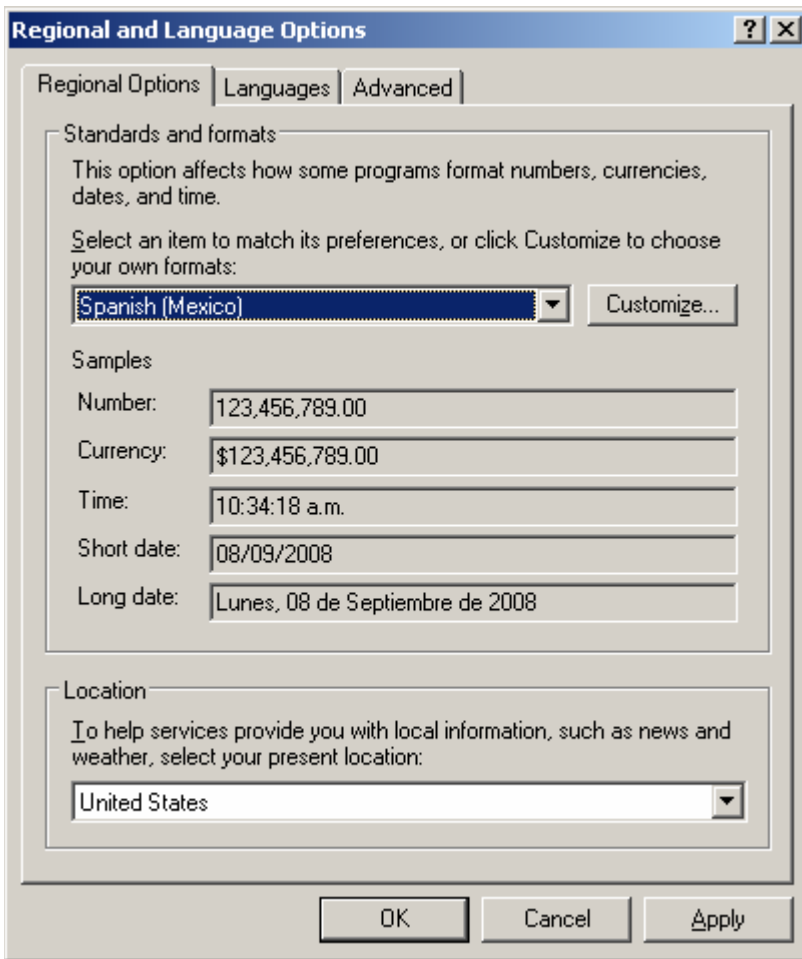

5. The “Language Bar” appears in the right side of the Taskbar to the left of the “System Tray,” displaying the language selection (in this example, EN = English - United States).




6. Click on the language selection to change it (in this example, ES = Spanish - Mexico).




7. Begin the application installation.




8. When finished, return to the control panel and select “Regional and Language Options.”

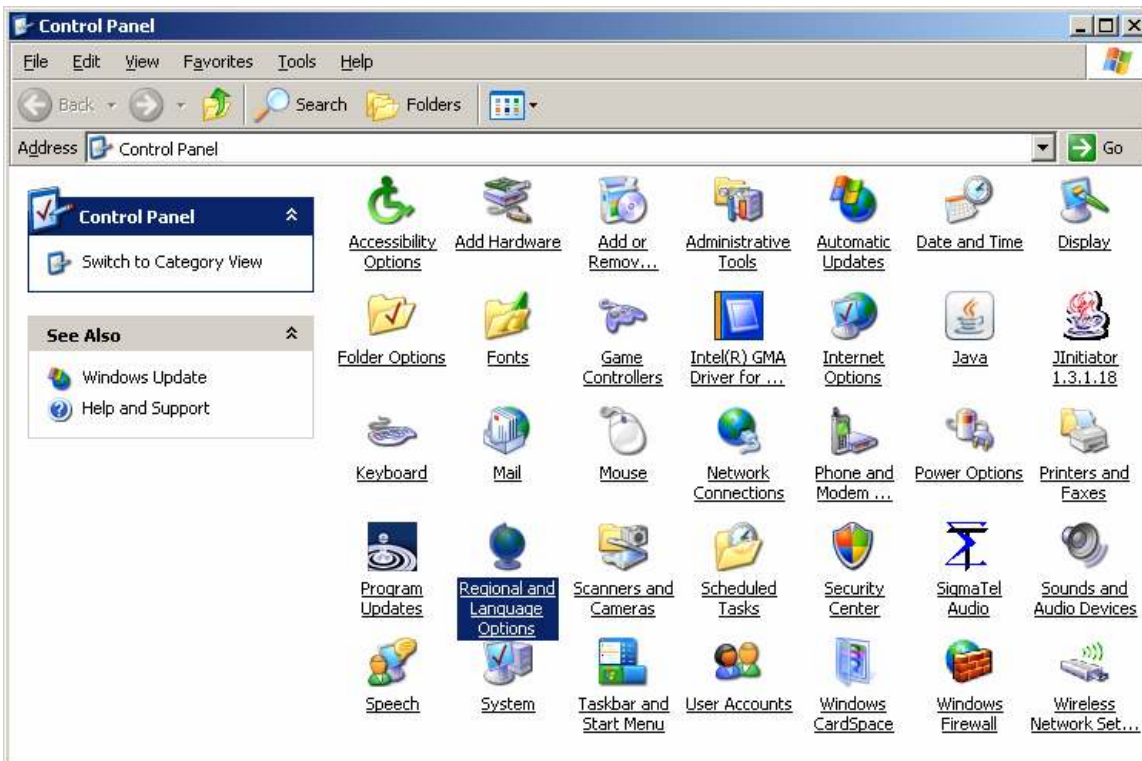


9. Click on the “Regional Options” tab and change the language selection back to its former value.

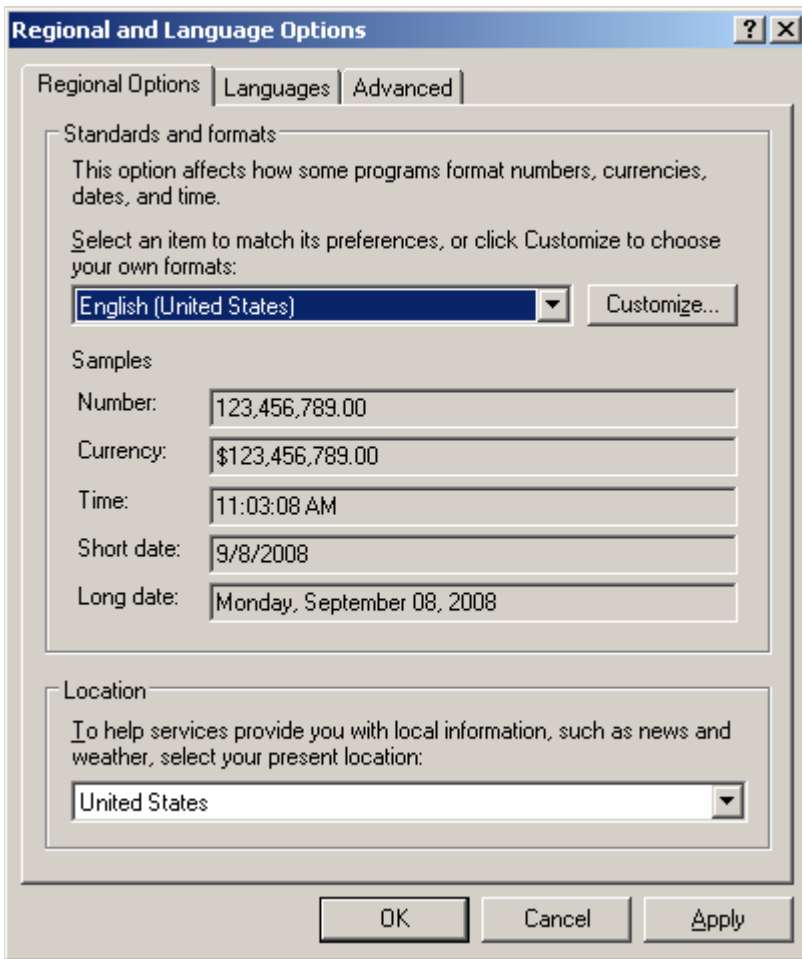


10. Click on the “Languages” tab and then click on the “Details” button.

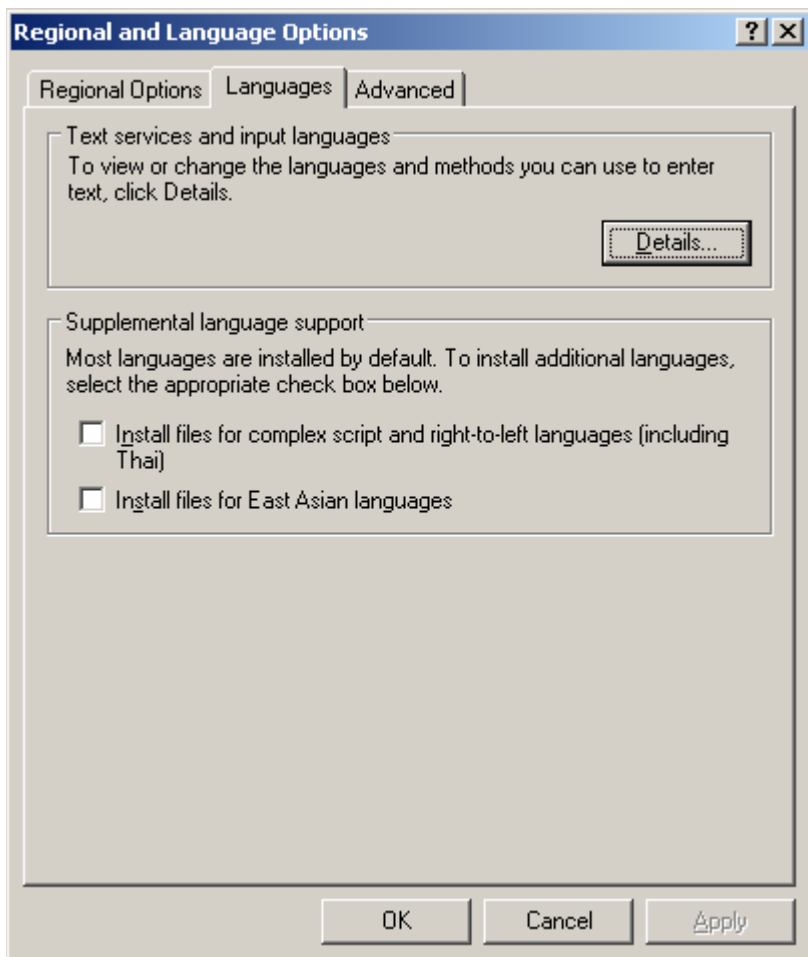

11. Click on the “Language Bar” button near the bottom of the panel.

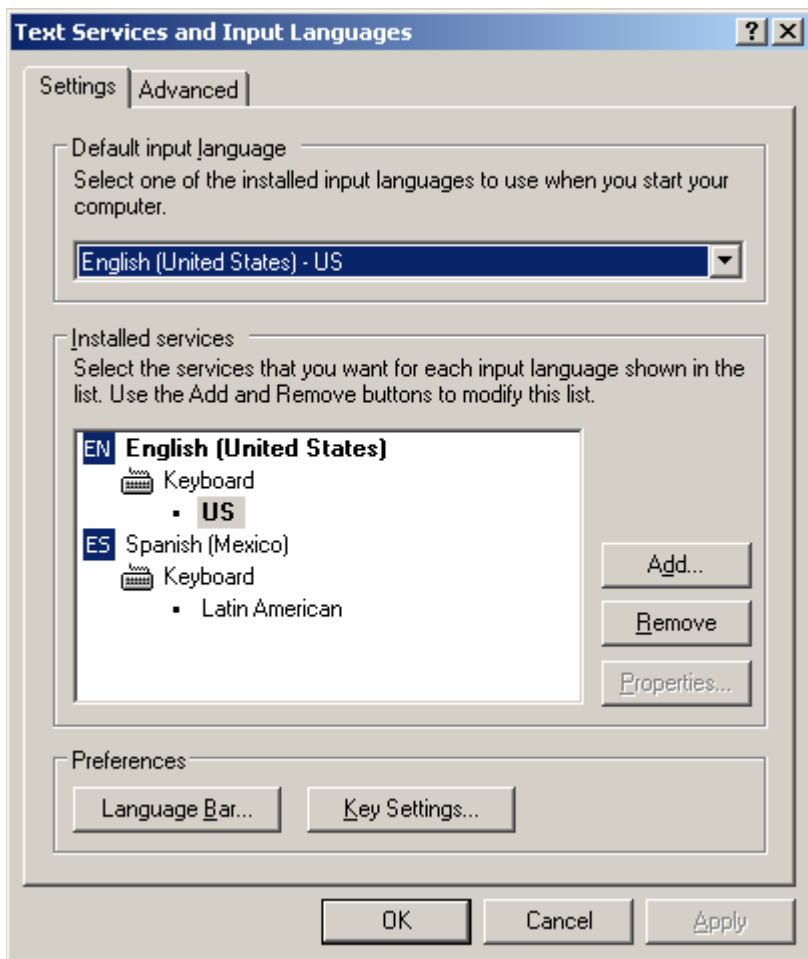

12. Uncheck the checkbox that reads “Show the Language bar on the desktop.” This will remove the language bar from the Taskbar.

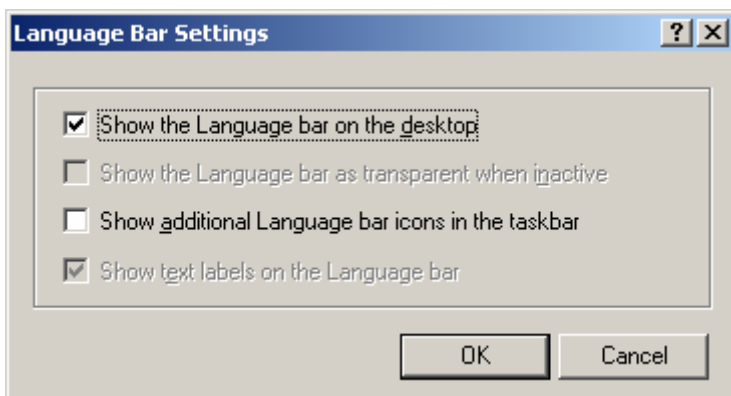

13. Click on the “OK” button and then close the “Control Panel.”

